



EZ BioPac

8 ADVANTAGES, ONE SMART SOLUTION

- EASE OF FILL:** Traditional single port powder bags have narrow openings, slowing fill time. EZ BioPac's generous opening speeds filling while reducing risk of overfills, spills or surface contamination.
- FINE-TUNING:** Narrow-neck bags make removing material to fine-tune weight difficult. EZ BioPac's big opening allows fast fine-tuning with a large scoop.
- 71% FASTER FILLS:** In direct-comparison trials, the EZ BioPac bag's larger opening and skirt reduced fill times substantially, compared to traditional 2D transfer bags.

FILLING TRIALS: Time to Fill to 5.0 kg		EZ BIOPAC®	2D TRANSFER BAG
TEST #	TEST MEDIA	PROCESS FILL TIME (MIN:SEC)	PROCESS FILL TIME (MIN:SEC)
Trial #1	Magnesium Sulfate	2:03.91	3:09.21
Trial #2	Magnesium Sulfate	1:54.66	3:01.06
Trial #3	Magnesium Sulfate	1:37.07	3:24.13

- REDUCED CONTAMINATION RISK:** Competitive single-use bag surfaces can become contaminated from product spills or overflows. EZ BioPac's outer skirt folds down to protect the frame and bag exterior, then up again to seal contaminants inside.
- CONVENIENT SEALING OPTIONS:** EZ BioPac bags can be sealed securing with attached cable ties.
- SEPARATE DISCHARGE OUTLET:** Unlike competitive 2D bags, EZ BioPac bags feature a separate discharge outlet that ensures a clear, clean discharge path and fast product transfer.
- 18% FASTER DISCHARGE:**

DISCHARGE TRIALS — Time to Discharge 5.0 kg		EZ BIOPAC®	2D TRANSFER BAG
TEST #	TEST MEDIA	DISCHARGE TIME (MIN:SEC)	DISCHARGE TIME (MIN:SEC)
Trial #1	Flour	1:49.0	1:45.0
Trial #2	Flour	1:51.81	2:03.07
Trial #3	Flour	1:21.42	2:10.70

- NO PRODUCT RESIDUE:** EZ BioPac's ArmorFlex®114 film has static dissipative properties, preventing product from clinging inside the bag, maximizing value and ensuring accurate dispense volumes.

5 Liters

Tri-Clamp Size	Tri-Clamp Type	Neck Style	ISO 7 Build	Gamma Irradiated
3"	Non-Gasket	V-Fold	EZ005V3A	EZ005V3A-G
3"		Weloc	EZ005W3A	EZ005W3A-G
4"	Non-Gasket	V-Fold	EZ005V4A	EZ005V4A-G
4"		Weloc	EZ005W4A	EZ005W4A-G

10 Liters

Tri-Clamp Size	Tri-Clamp Type	Neck Style	ISO 7 Build	Gamma Irradiated
3"	Non-Gasket	V-Fold	EZ010V3A	EZ010V3A-G
3"		Weloc	EZ010W3A	EZ010W3A-G
4"	Non-Gasket	V-Fold	EZ010V4A	EZ010V4A-G
4"		Weloc	EZ010W4A	EZ010W4A-G

25 Liters

Tri-Clamp Size	Tri-Clamp Type	Neck Style	ISO 7 Build	Gamma Irradiated
3"	Non-Gasket	V-Fold	EZ025V3A	EZ025V3A-G
3"		Weloc	EZ025W3A	EZ025W3A-G
4"	Non-Gasket	V-Fold	EZ025V4A	EZ025V4A-G
4"		Weloc	EZ025W4A	EZ025W4A-G

50 Liters

Tri-Clamp Size	Tri-Clamp Type	Neck Style	ISO 7 Build	Gamma Irradiated
4"	Non-Gasket	V-Fold	EZ050V4A	EZ050V4A-G
4"		Weloc	EZ050W4A	EZ050W4A-G
6"	Sanitary	V-Fold	EZ050V6B	EZ050V6B-G
6"		Weloc	EZ050W6B	

100 Liters

Tri-Clamp Size	Tri-Clamp Type	Neck Style	ISO 7 Build	Gamma Irradiated
6"	Sanitary	V-Fold	EZ100V6B	EZ100V6B-G
6"		Weloc	EZ100W6B	

ACCESSORIES

Fill Frames

Description	Part Number
Frame, EZ BioPac, 5/10L	HW10429-01
Frame, EZ BioPac, 25/50L	HW10376-02
Frame, EZ BioPac, 100L	HW10428-01
Frame, EZ BioPac, 1L, HDPE/SS	HW12137-01

Lifting Bars

Description	Part Number
Lifting Bar, 4 Loop, 14" Dia Body	HW11200-01
Lifting Bar, 4 Loop, 14" Dia Body, Low Profile	HW11200-02
Lifting Bar, 8 Loop, 20" Dia Body	HW11089-01
Lifting Bar, 8 Loop, 20" Dia Body, Low Profile	HW11089-02
Hook, EZ BioPac, 5L/10L	HW12017-01



SCAN ME



BEYOND BOUNDARIES™

ILC Dover LP
 One Moonwalker Road
 Frederica, DE 19946 USA
 +1.302.335.3911
 +1.800.631.9567
 customer_service@ilcdover.com

www.ilcdover.com

